

Спектрометры портативные Рамановские серии ATR3110PS

Технические характеристики

По вопросам продаж и поддержки обращайтесь:

Алматы (727)345-47-04
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Саранск (8342)22-96-24
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(727)345-47-04

Беларусь +375-257-127-884

Узбекистан +998(71)205-18-59

Киргизия +996(312)96-26-47

эл.почта: otp@nt-rt.ru || сайт: <https://optosky.nt-rt.ru/>

Datasheet

High-sensitivity High-resolution Portable Raman Spectrometer

ATR3110PS

Feature:

- Ultra-high sensitivity FFT-CCD TE-cooled;
- low noise circuit;
- Powerful embedded software;
- Fluorescent background eliminate;
- Peak finding and display;
Win 10 operation system;
- USB 2.0;
- User friendly human-machine interface;
- Remote control via LAN;

Application:

- Biological science
- Pharmaceutical engineering
- Forensic analysis
- Agriculture and food safety
- Gemstone
- Environmental science

Description:

ATR3110PS Raman Spectrometer is TE-cooled, high-sensitivity, enhanced designed for broadband ranges. all of optical path, PCB, signal processing method have been made optimized processed to obtain >15times higher SNR than ATR20007, nearly 100 times higher than 2000cm⁻¹. ATR3110PS employs low noise CCD signal process circuit, noise<3counts.

ATR3110PS employs 110/220V power supply, DC supply via 5V adaptor. Easy to take and field operation.

PN	Wavelength (nm)	Wavenumber range cm ⁻¹
ATR3110PS-473	473	200-4000
ATR3110PS-532	532	200-3600
ATR3110PS-785-27	785	200-2700
ATR3110PS-785-40		200-4300
ATR3110PS-830	830	200-4000
ATR3110PS-1064	1064	200-2600
Available in custom made wavelength		

Remark:

- Measuring method is based on ASTM E2529-06;
- Available in custom design, resolution can be increased by around 1/3, resulting in lower sensitivity;



1 Specifications

ATR3110PS System			
Operating system	Windows		
Integration time	6ms-30mins		
Power voltage	DC 5V(+/-5%)		
Operating Temp	-10~40 °C		
Operating humidity	< 95%		
Dimension(L*W*H)	30×22.5×13.2 cm		
Weight	5.5 Kg		
Reliability			
Spectral stability	$\sigma/\mu < 0.5\%$ (COT 8 hours)		
Temp stability	Spectral shift $\leq 1 \text{ cm}^{-1}$ (10-40 °C)		
Variation of intensity (in 5 ~ 40 °C)	<±5%		
Optical parameters			
Spectral range (cm ⁻¹)	200-2700	200-3500	200-4300
resolution (cm ⁻¹) / 50μm slit size	6.5nm	9nm	12nm
SNR	>3000:1 (918 cm ⁻¹ of Acetonitrile, 10s accumulation, 200mW)		
Sensitivity	3200:1		
Optical system	f/4 C-T crossed optical path		
focusing	98 mm for incidence and output		
Detector			
Item	Ultra-high sensitivity, quick cooling CCD		
Detector cooled down to	-5 °C		
Detecting range	200-1100 nm		
Effective pixels	2048*64		
Dynamic range	50000: 1		
Pixel size	14μm×200μm		
Exciting Laser			
Central wavelength	785nm (+/-1nm)		
FWHM	0.08 nm		
Power output	≥500 mW		
Power stability	$\sigma/\mu < \pm 0.2\%$		
Raman probe			

Datasheet

Operating distance	6 mm
Rayleigh scattering resistance	OD>8
Numerical Aperture	0.3
Aperture	7mm

2 Optical Performance

2.1 General spectral performance

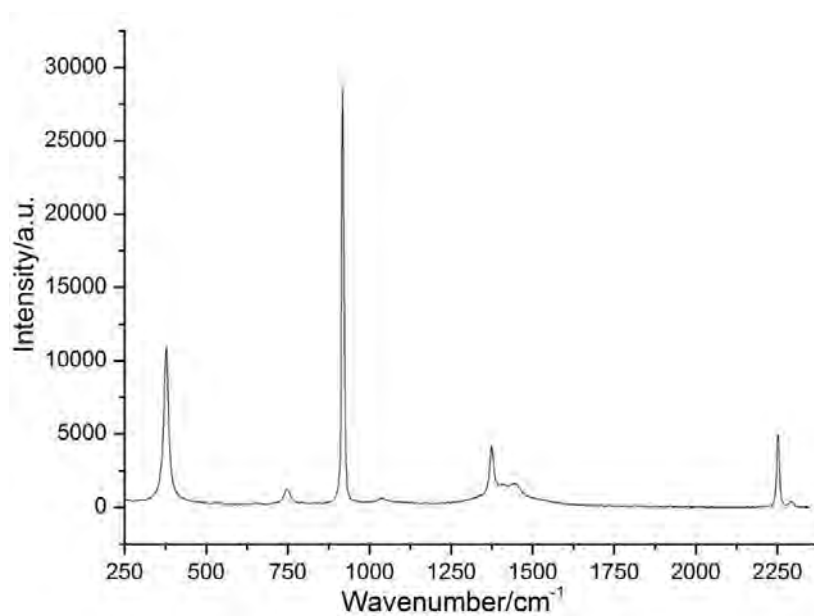


Figure 1 Raman spectra of acetonitrile

Datasheet

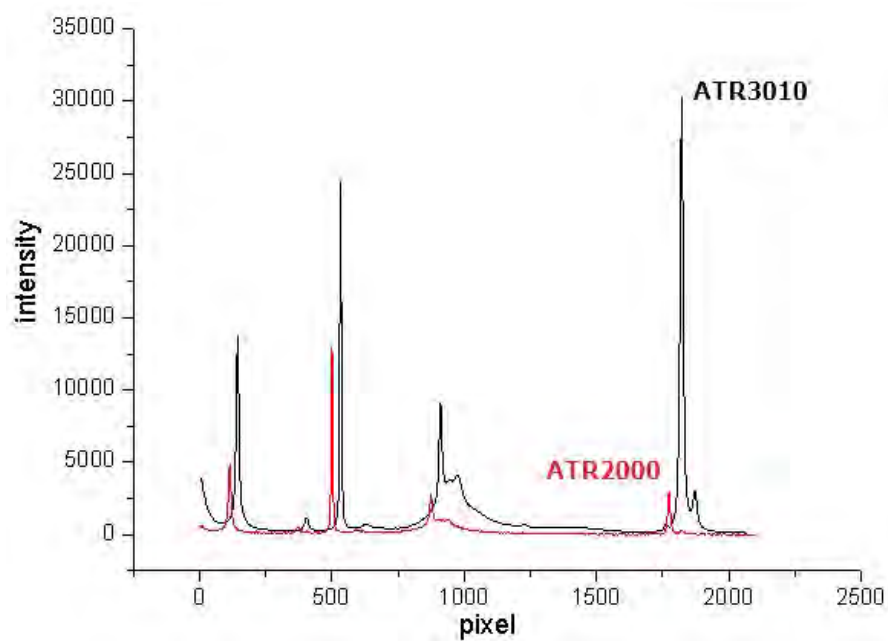


Figure 2 Sensitive of ATR3000 vs ATR2000

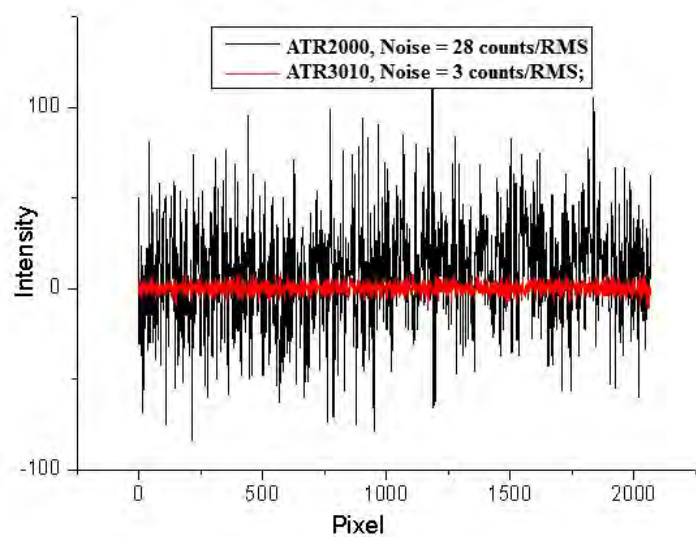


Figure 3 Noise of ATR3000 vs ATR2000

2.2 Spectral Resolution

2.2.1 Raman spectral of Tylenol

Laser Power 200 mW
Integration time: 10 s
Filter level 1

Raman spectra of Tylenol showed the resolution condition in the long wavelength region. That is better than 6 cm^{-1} .

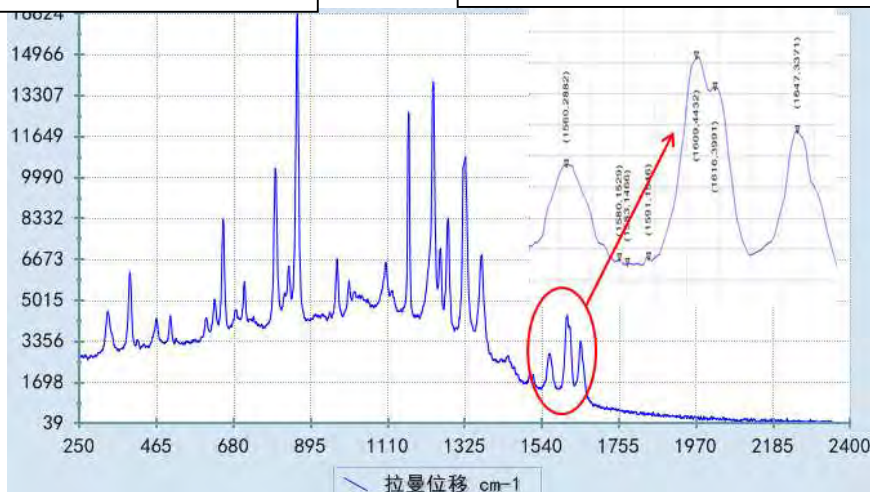


Fig.2.2 Raman spectrum of Tylenol, the vibration mode $1610/1615\text{ cm}^{-1}$ can be resolved.

2.2.2 Raman spectral of petrol

Laser Power 200 mW
Integration time: 10 s
Filter level 1

Raman spectra of petrol 93# showed the resolution condition in the short wavelength region.

Datasheet


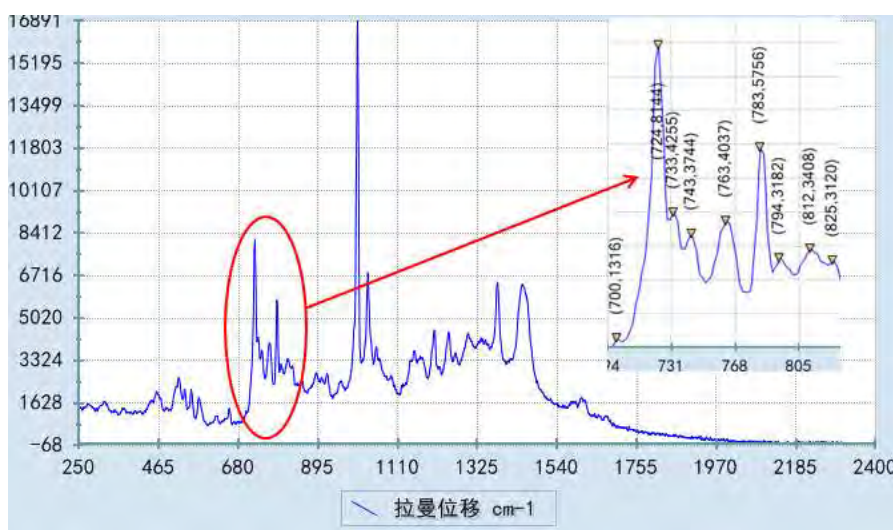


Fig.2.3 Raman spectrum of petrol 93#, the vibration mode 723/732/742 cm^{-1} can be resolved.

3 Reliability

Figure 3.1 and Figure 3.2 showed the temperature reliability testing results of five ATR3000 portable Raman spectrometers. The testing temperature range was from 5 °C to 40 °C. The spectrometer was kept more than 1 hour at every temperature spots. Acetonitrile was used as the standard sample in the testing. The testing results were calculated using 918 cm^{-1} of acetonitrile. The wavenumber shift was 1 cm^{-1} or less (as show in Fig. 3.1). The peak intensity variation was less than 10% (as show in Fig. 4).

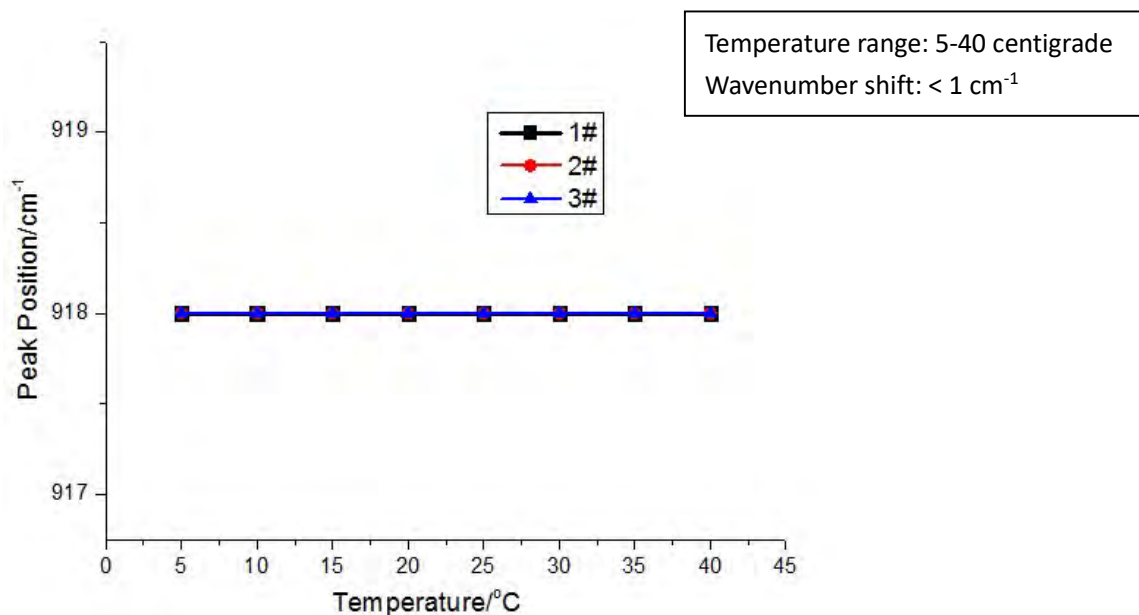

Datasheet

Fig. 3.1 Wavenumber shift results testing from 5 °C to 40 °C of fives ATR3000 portable Raman spectrometers

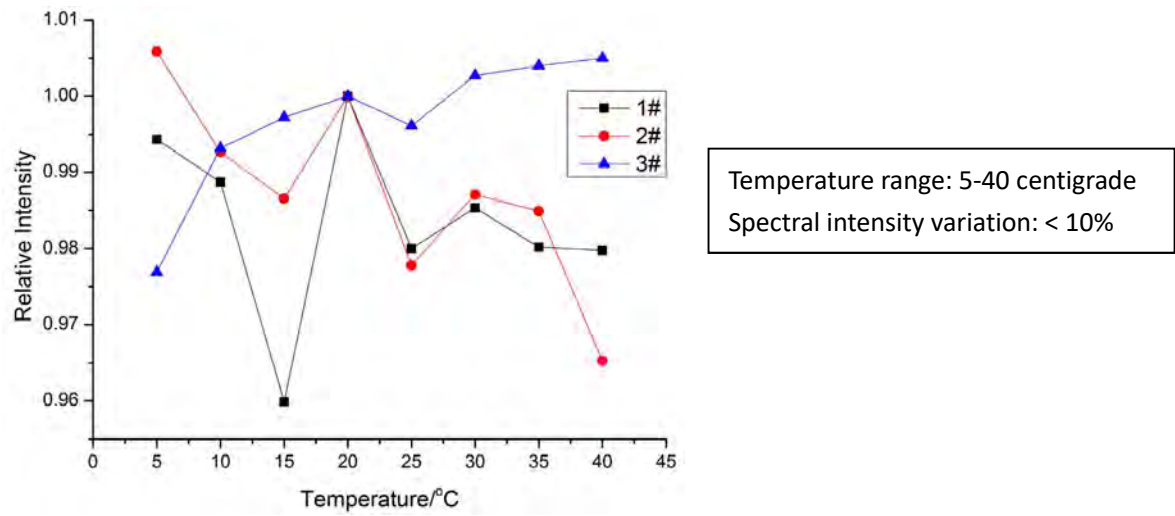


Figure 4 Intensity variation testing from 5 °C to 40 °C of three ATR3000 portable Raman spectrometers

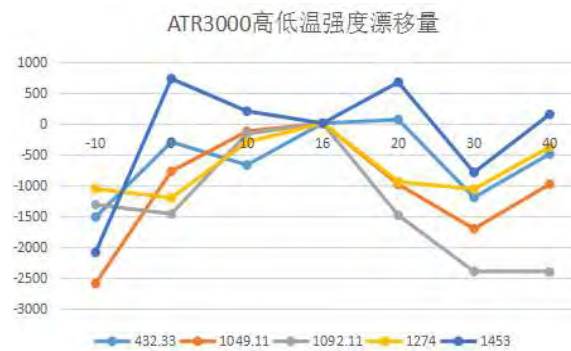


Figure 5 Intensity variation -10 °C to 40 °C of ATR3000 portable Raman spectrometers, sample is alcohol.

Datasheet

2. Measuring accessories



Fig 2 Solid, powder measuring probe



Fig 1 Fluid sample cell (Thermo bottle)



Fig 2 Fluid sample cell (Liquid chromatography bottle) **(Optional)**

Datasheet



Fig 5 Raman probe gun (optional)



Fig 6 Measuring adjustable holder (Optional)

3. Other excitation wavelength:

Datasheet

ITEM No.	Excitation Wavelength (nm)	Maximum laser power (mW)	Spectral range (cm ⁻¹)	resolution (cm ⁻¹)	Feature
ATRPS31100-27	785	550	200-2700	6	Available for most application
ATRPS31100-35			200-3500	8	
ATRPS31100-43			200-4300	10	
ATR3110PS-1064	1064	500	200-2600	13	Fluorescence-free, non-destructive ,high-sensitivity, high-SNR, Available samples: dark-color samples, fluorescence sample, biology sample, bacteria, plastic, fuel, petroleum product, vegetable oil, explosive etc.
ATR3110PS-830	830	550	200-3300	7	higher skin permeance suit to biological samples, eg. Non-invasive blood glucose, early cancer diagnosis
ATR3110PS-266	266	50	200-3000	25	
ATR3110PS-532	532	100	200-3600	10	
ATR3110PS-633	633	80	200-3200	10	



По вопросам продаж и поддержки обращайтесь:

Алматы (727)345-47-04
Ангарск (3955)60-70-56
Архангельск (8182)63-90-72
Астрахань (8512)99-46-04
Барнаул (3852)73-04-60
Белгород (4722)40-23-64
Благовещенск (4162)22-76-07
Брянск (4832)59-03-52
Владивосток (423)249-28-31
Владикавказ (8672)28-90-48
Владимир (4922)49-43-18
Волгоград (844)278-03-48
Вологда (8172)26-41-59
Воронеж (473)204-51-73
Екатеринбург (343)384-55-89

Иваново (4932)77-34-06
Ижевск (3412)26-03-58
Иркутск (395)279-98-46
Казань (843)206-01-48
Калининград (4012)72-03-81
Калуга (4842)92-23-67
Кемерово (3842)65-04-62
Киров (8332)68-02-04
Коломна (4966)23-41-49
Кострома (4942)77-07-48
Краснодар (861)203-40-90
Красноярск (391)204-63-61
Курск (4712)77-13-04
Курган (3522)50-90-47
Липецк (4742)52-20-81

Магнитогорск (3519)55-03-13
Москва (495)268-04-70
Мурманск (8152)59-64-93
Набережные Челны (8552)20-53-41
Новокузнецк (3843)20-46-81
Ноябрьск (3496)41-32-12
Новосибирск (383)227-86-73
Омск (3812)21-46-40
Орел (4862)44-53-42
Оренбург (3532)37-68-04
Пенза (8412)22-31-16
Петрозаводск (8142)55-98-37
Псков (8112)59-10-37
Пермь (342)205-81-47

Ростов-на-Дону (863)308-18-15
Рязань (4912)46-61-64
Самара (846)206-03-16
Санкт-Петербург (812)309-46-40
Саратов (845)249-38-78
Севастополь (8692)22-31-93
Саранск (8342)22-96-24
Симферополь (3652)67-13-56
Смоленск (4812)29-41-54
Сочи (862)225-72-31
Ставрополь (8652)20-65-13
Сургут (3462)77-98-35
Сыктывкар (8212)25-95-17
Тамбов (4752)50-40-97
Тверь (4822)63-31-35

Тольятти (8482)63-91-07
Томск (3822)98-41-53
Тула (4872)33-79-87
Тюмень (3452)66-21-18
Улан-Удэ (3012)59-97-51
Уфа (347)229-48-12
Хабаровск (4212)92-98-04
Чебоксары (8352)28-53-07
Челябинск (351)202-03-61
Череповец (8202)49-02-64
Чита (3022)38-34-83
Якутск (4112)23-90-97
Ярославль (4852)69-52-93

Россия +7(495)268-04-70

Казахстан +7(727)345-47-04

Беларусь +375-257-127-884

Узбекистан +998(71)205-18-59

Киргизия +996(312)96-26-47

эл.почта: otp@nt-rt.ru || сайт: <https://optosky.nt-rt.ru/>